



COLOR TOUCH SCREEN DISPLAY

SAFE HEAT MODE TO PROTECT AND  
EXTEND FILAMENT LIFE


PIPETTE COOKBOOK PROGRAM DIRECTORY

LINE REPEAT MODE SIMPLIFIES  
MULTI-LINE PROGRAMMING

HELP TOPICS AND ERROR DETECTION

GLOSSARY WITH MICROPIPETTE AND  
PULLER TECHNOLOGY

RECORD OF LAST TWO PULL RESULTS

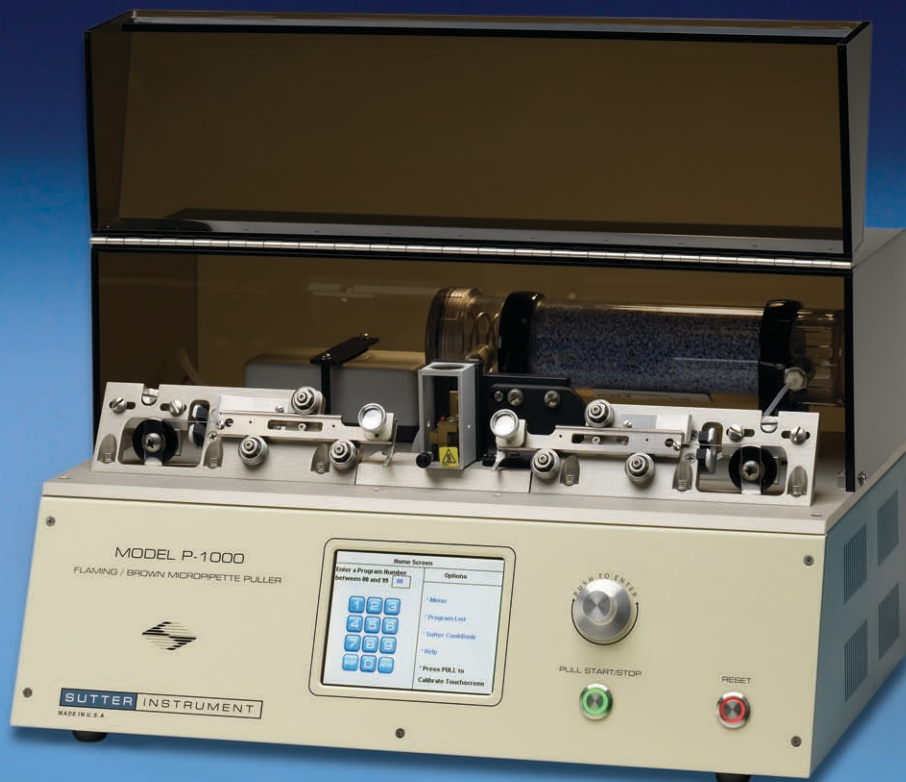

## P-1000 NEXT GENERATION MICROPIPETTE PULLER

Developed through years of experience with the Flaming-Brown style micropipette pullers, and infused with leading-edge technology, Sutter is proud to introduce the **P-1000** micropipette puller. The most obvious new feature is the color touch-screen display that provides an intuitive and full featured interface.

The extensive library of programs found in the popular Sutter Pipette Cookbook has been incorporated into the **P-1000** puller and are available to the user. You need only specify the glass, filament, and type of pipette you require, and a suitable program is identified and available for installation. This takes the guesswork out of pipette pulling and simplifies programming.

Copy and Paste functions assure that programs can be easily written and the line repeat mode simplifies multi-line programming. A safe heat mode is an additional feature that helps the user avoid using

heat settings that might damage or burn out the filament. When the safe heat mode is turned on, the puller will "check" the installed heat and alert the user when a given heat value is too low or too high in relation to the ramp value. User notes can be added to each program for annotating important information.

New features for the **P-1000** include: diagnostic testing of all puller components, built-in error detection of air pressure loss or filament burnout, easy access to ramp test, measurement of jaw temperature, and access to previous pull results with the heat-on times for each cycle of the program. Help topics are preloaded to assist with on-site troubleshooting, and the built-in glossary includes text, pictures, and diagrams explaining the terms used in micropipette fabrication. A rotary dial is offered as an alternative to the touch pad for numerical entry.

### **P-1000**

Flaming/Brown type micropipette puller, manual

Each puller comes with an assortment of 4 filaments and a sample box of BF100-50-10 and BF150-86-10 glass. Sutter pre-programs the P-1000 with a 2.5mm box filament unless an alternative filament is requested.

**SUTTER INSTRUMENT**

One Digital Drive • Novato • CA 94949 • Phone 415.883.0128  
Fax 415.883.0572 • Web [www.sutter.com](http://www.sutter.com) • Email [info@sutter.com](mailto:info@sutter.com)